



SERVICE REPORT

 MIG-315, Balaji Nagar, Kukatpally, Hyderabad - 500 072, Ph (O) : 9505555984, E-mail : krishna@tecpharma.in	Date : 24/11/2020	S/R : 1610
	Engineer's Name : D. Ashole	
	Mobile No. : 8897885376	
Customer's Name :	Type of Services	
Address Indian Immunological Limited	1. Installation <input checked="" type="checkbox"/>	6. Paid Service <input type="checkbox"/>
Rakshapuram, Gachibowli post	2. Breakdown <input type="checkbox"/>	7. Calibration <input type="checkbox"/>
Hyderabad-500032	3. AMC <input type="checkbox"/>	8. IQ/OQ/PQ <input checked="" type="checkbox"/>
Telangana, India	4. Warranty <input type="checkbox"/>	9. Application <input type="checkbox"/>
	5. Inspection <input type="checkbox"/>	10. Goodwill <input type="checkbox"/>
P.O. No.	Job Completed : Yes <input checked="" type="checkbox"/> No <input type="checkbox"/>	
E-mail :		
Instrument	Model No.	Sr. No. / INST ID
Haier Ultra Low Temperature Freezer	DW-86L729BP	BEGY8EBA009GL880001/ B5/ULF/218
Problem As Reported by the Customer : Haier ULT Freezer Installation.		
Service Done : Haier Ultra Low Temperature freezer installation completed, performed IQ, OQ and PQ Qualifications. During 72 hours temperature mapping without load, 72 hours temperature mapping with load at $-78 \pm 1^{\circ}\text{C}$ power failure study and door open study done. now ultra low temperature		
Note : freezer working satisfactorily.		
Spare Parts Used / Required	To Be filled by Customer	
NA	<input type="checkbox"/> Instrument Installed & Found Working Satisfactorily <input type="checkbox"/> Prevention Maintenance Carried Out Satisfactorily <input type="checkbox"/> Fault Rectified & Instrument Working Satisfactorily <input type="checkbox"/> Please await our written approval for parts replacement <input type="checkbox"/> Calibration done Satisfactorily <input type="checkbox"/> Comments (If Any)	
Engineer's Signature : D. Ashole 24/11/2020	Name : P. Sukumar	[Signature] Customer Signature
	Date : 24-11-20	* Only for parts Changed & all other charges will be extra as applicable



Note : Parts replaced are chargeable except during warranty period. Consumables used for the instrument are not covered by warranty & hence are chargeable. Parts replaced due to mishandling during operation or otherwise will also be charged in every case. In case of any difficulty or suggestion, please contact Service Manager.

SERVICE REPORT

 MIG-315, Balaji Nagar, Kukatpally, Hyderabad - 500 072, Ph (O) : 9505555984, E-mail : krishna@tecpharma.in	Date : 24/11/2020.	S/R : 1611
	Engineer's Name : D. Ashok	
	Mobile No. : 8897885376	
Customer's Name :	Type of Services	
Address Indian Immunological Limited	1. Installation <input checked="" type="checkbox"/>	6. Paid Service <input type="checkbox"/>
Raothapuram, Gachibowli post	2. Breakdown <input type="checkbox"/>	7. Calibration <input type="checkbox"/>
Hyderabad - 500032	3. AMC <input type="checkbox"/>	8. IQ/OQ/PQ <input checked="" type="checkbox"/>
Telangana, India	4. Warranty <input type="checkbox"/>	9. Application <input type="checkbox"/>
	5. Inspection <input type="checkbox"/>	10. Goodwill <input type="checkbox"/>
P.O. No.	Job Completed : Yes <input checked="" type="checkbox"/> No <input type="checkbox"/>	
E-mail :		
Instrument	Model No.	Sr. No. / INST ID
Haier Ultra Low Temperature Freezer	DW-86L729BP	BEO6Y8EBA00Q6L880006/ B5/ULF/219
Problem As Reported by the Customer : Haier ULT Freezer installation		
Service Done : Haier Ultra Low Temperature Freezer installation, completed, performed IQ, OQ and PQ qualification, during 72 hours temperature mapping without load, 72 hours temperature mapping with load at $-70^{\circ}\text{C} \pm 1^{\circ}\text{C}$, power failure study and Door open study done, now		
Note : the Ultra Low temperature freezer working satisfactorily.		
Spare Parts Used / Required	To Be filled by Customer	
/	<input type="checkbox"/> Instrument Installed & Found Working Satisfactorily <input type="checkbox"/> Prevention Maintenance Carried Out Satisfactorily <input type="checkbox"/> Fault Rectified & Instrument Working Satisfactorily <input type="checkbox"/> Please await our written approval for parts replacement <input type="checkbox"/> Calibration done Satisfactorily <input type="checkbox"/> Comments (If Any)	
NA	Name : P. Subramaniam Date : 24-11-20	
/	Customer Signature * Only for parts Changed & all other charges will be extra as applicable	
Engineer's Signature : D. Ashok 24/11/2020		

Note : Parts replaced are chargeable except during warranty period. Consumables used for the instrument are not covered by warranty & hence are chargeable. Parts replaced due to mishandling during operation or otherwise will also be charged in every case. In case of any difficulty or suggestion, please contact Service Manager.

SERVICE REPORT

 MIG-315, Balaji Nagar, Kukatpally, Hyderabad - 500 072, Ph (O) : 9505555984, E-mail : krishna@tecpharma.in	Date : 24/11/2020	S/R : 1612
	Engineer's Name : D. Ashok	
	Mobile No. : 8897885376	
Customer's Name :	Type of Services	
Address Indian Immunological Limited	1. Installation <input checked="" type="checkbox"/>	6. Paid Service <input type="checkbox"/>
Rakshapuram, Gachibowli post	2. Breakdown <input type="checkbox"/>	7. Calibration <input type="checkbox"/>
Hyderabad - 500032	3. AMC <input type="checkbox"/>	8. IQ/OQ/PQ <input checked="" type="checkbox"/>
Telangana, India	4. Warranty <input type="checkbox"/>	9. Application <input type="checkbox"/>
	5. Inspection <input type="checkbox"/>	10. Goodwill <input type="checkbox"/>
P.O. No.	Job Completed : Yes <input checked="" type="checkbox"/> No <input type="checkbox"/>	
E-mail :		
Instrument	Model No.	Sr. No. / INST ID
Haier Ultra Low temperature Freezer	DW-P6L729BP	BEOGY8EBA00QGLP0007/ B5/ULF/220
Problem As Reported by the Customer : Haier ULT Freezer installation		
Service Done : Haier Ultra Low Temperature Freezer Installation, Completed, performed IQ, OQ and PQ Qualification during 72 hours temperature mapping without load, 72 hours temperature mapping with load at $-70^{\circ}\text{C} \pm 1^{\circ}\text{C}$, power failure study and Door open study done.		
Note : now the Ultra Low temperature freezer working satisfactorily		
Spare Parts Used / Required	To Be filled by Customer	
NA	<input type="checkbox"/> Instrument Installed & Found Working Satisfactorily <input type="checkbox"/> Prevention Maintenance Carried Out Satisfactorily <input type="checkbox"/> Fault Rectified & Instrument Working Satisfactorily <input type="checkbox"/> Please await our written approval for parts replacement <input type="checkbox"/> Calibration done Satisfactorily <input type="checkbox"/> Comments (If Any)	
Engineer's Signature : D. Ashok 24/11/2020	Name : P. Sukuman Date : 24.11.20	 Customer Signature * Only for parts Changed & all other charges will be extra as applicable

Note : Parts replaced are chargeable except during warranty period. Consumables used for the instrument are not covered by warranty & hence are chargeable. Parts replaced due to mishandling during operation or otherwise will also be charged in every case. In case of any difficulty or suggestion, please contact Service Manager.